

Operating Instructions

Sartobind STIC[®] PA

Void Volume Optimized Capsules and Cassettes
With 4 mm Bed Height



4 mm: nano
1 mL

75 mL

200 mL

400 mL

600 mL

Jumbo 2.5 L


Cassette 0.8 L


1000032238




SARTORIUS


Read operational instructions carefully before using Sartobind capsules.

 Use of the products in applications not specified or not described in this manual, may result in improper function, personal injury, or damage of the product or material. The products are supplied as non-sterile unless otherwise expressly described. The membrane is dried from glycerol.

 Die Verwendung dieser Produkte für Anwendungen, für die sie nicht bestimmt oder nicht in dieser Anleitung beschrieben sind, können zu einer schlechteren Funktion, Zerstörung der Produkte oder sogar zu Verletzungen von Mensch und Material führen. Die Produkte sind nicht steril sofern dies nicht ausdrücklich anders beschrieben ist. Die enthaltene Membran wird aus Glycerin getrocknet.

- ⚠ L'utilisation des produits pour des applications non-spécifiées ou décrites dans ce manuel peut causer un dysfonctionnement, une destruction du produit, des dommages matériels ou même corporels. Les produits sont fournis non-stériles, sauf indication contraire expressément mentionnée. La membrane est séchée avec de la Glycérine.
- ⚠ La utilización de este producto en aplicaciones ajenas o no establecidas en el manual de operación, puede provocar un mal funcionamiento del producto, del material, así como daños personales. Los productos suministrados no son estériles a menos que se describa lo contrario. La membrana ha sido secada de glicerina.

 把产品用于手册中没指定或描述的应用, 可能导致产品失效, 人员受伤, 或使产品及物料受损。囊式膜柱以非灭菌的形式供应。膜用甘油脱水。

 当製品を該当しない用途、あるいは当製品取扱説明書に記載されていない応用分野において使用した場合、当製品の機能上の不具合や損傷、人体への危害、あるいは他の物品の損傷を招く恐れがあります。特に明記のない場合、当製品は滅菌処理されていません。当メンブレンはグリセリンを用いて乾燥させてあります。

Intended use

The membrane chromatography products – also described as membrane adsorbers – should be used only once for flowthrough (negative) chromatography applications to avoid carryover as well as tedious and costly cleaning validation procedure.

Sartobind STIC® PA nano 1 mL has been developed as a scouting device for working with small sample volumes while retaining the cylindrical design of the large scale Membrane Adsorbers.

Sartobind STIC® PA 75 mL up to Jumbo 2.5 L have been developed for intermediate and pilot scale up to production scale in the biopharmaceutical industry.

Sartobind 0.8 L cassettes are used in the Pilot Scale Filter Holder of up to 10.4 L membrane volume for the biopharmaceutical production.

Table of Contents

1	Storage conditions	10
2	Introduction	11
3	Technical Data	16
4	Materials	18
5	Binding capacity	19
6	Installation	20
7	Operation	24
7.1	Venting	24
7.2	Cleaning and equilibration	27
7.3	Autoclaving	27
7.4	Recommended flow rates and equilibration volumes	28
7.5	Buffer conditions	30
7.6	Selection of pH and salt conditions	31
7.7	Contaminant removal from therapeutic proteins and other sources in flow-through mode	32

7.8	Sample preparation.....	33
7.9	Washing.....	33
7.10	Elution.....	34
7.11	Draining.....	34
7.12	Regeneration and Storage.....	35
7.13	Chemical stability.....	35
7.14	Operation of the Sartobind nano with peristaltic pumps or liquid chromatography (LC) systems.....	35
7.15	Scaling up.....	36
8	Integrity test by diffusion.....	40
8.1	Installation for test.....	40
8.2	Operation of test.....	41
9	Troubleshooting.....	44
10	Quality assurance.....	49

11	Ordering information	50
11.1	Products.....	50
11.2	Accessories.....	52
12	Dimensions and Connections	54

1 Storage Conditions

Sartobind capsules should be stored clean dry and away from direct sunlight in the closed bag and box at room temperature.

2 Introduction

The capsules and cassettes with 4 mm bed height are salt tolerant weak anion exchange chromatography devices based on macroporous membranes. They can be used for chromatographic separations in the downstream processing of viruses and proteins. The ion exchange ligands are coupled to a membrane which is fitted into a plastic housing for quick handling, making ion exchange purification nearly as easy as filtration.

The devices are constructed with optimized fluid channels. The capsules a central core and the cassettes a spacer element to minimize void volume. To set up and operate the Sartobind Jumbo (2.5 liter membrane volume) we recommend the Jumbo trolley (see chapter "11.2 Accessories", page 52). These products are intended for single use to avoid carryover as well as tedious and costly cleaning validation procedures (see also section "7.12 Regeneration and Storage", page 35). They are applied for contaminant removal from proteins and viral | virus like particles (VLP) vaccines in flow-through mode (negative chromatography) to bind DNA, residual protein, host cell proteins (HCP), endotoxins and viruses.

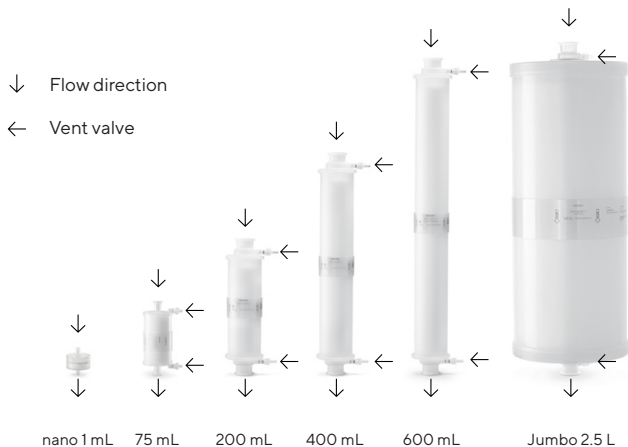


Fig. 1: Flow direction and position of vent valves of 4 mm capsules



Fig. 2: Flow direction and position of vent valve connection of 4 mm cassettes

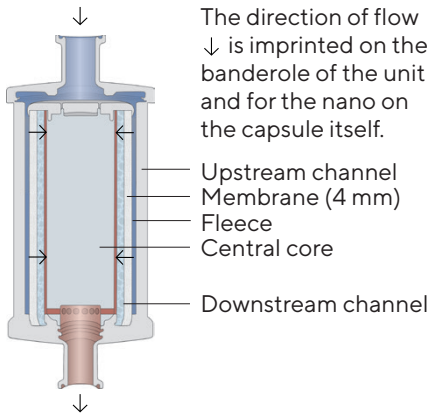



Fig. 3: Construction and flow path inside the capsules

 Devices should be visually inspected before use. In case of visible damage, the module must be replaced. Close vent valves of capsules before use by screwing the valve clockwise. For the cassettes close the clamps at the manifold set.

For the nano and 75 mL devices the central core is made from a solid polypropylene cylinder. For the larger capsules it is made from a self-contained air filled polypropylene cylinder. The interior of the core is inaccessible for gases and fluids. The two flat membrane stacks of the cassettes are separated by a central spacer element.



Fig. 4: Side view cassette; highlighted section see Fig. 5

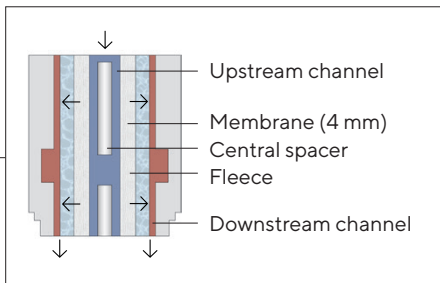


Fig. 5: Construction and flow path inside the 4 mm cassette

3 Technical Data

Membrane volume (MV)	1 mL	75 mL
Nominal membrane area	36.4 cm ²	2,700 cm ²
Bed height	4 mm	4 mm
Design	Cylindrical	Cylindrical
Sartobind STIC PA typical 10% dynamic binding capacity*	50 mg	3.75 g
Maximum pressure bar (MPa, psig) at 20°C	4 (0.4, 58)	4 (0.4, 58)
Maximum pressure during venting bar (MPa, psig) at 20°C	-	0.5 (0.05, 7)
Nominal void volume (mL)	3.5	200
Nominal void volume (MV)	3.5	2.4
Approximate weight	10 g	400 g

1 mL membrane = 36.4 cm² membrane

Ion capacity per cm² of membranes: 18–22 µeq

Short term pH stability 2–14 refers to cleaning in place procedure described in section “7.2 Cleaning and equilibration”, page 27

200 mL	400 mL	600 mL	2.5 L	0.8 L
7,300 cm ²	14,600 cm ²	22,000 cm ²	91,000 cm ²	29,000 cm ²
4 mm	4 mm	4 mm	4 mm	4 mm
Cylindrical	Cylindrical	Cylindrical	Cylindrical	Flat sheet
10 g	20 g	30 g	127 g	40
4 (0.4, 58)	4 (0.4, 58)	4 (0.4, 58)	3 (0.3, 43,5)	2 (0.2, 29)
0.5 (0.05, 7)	0.5 (0.05, 7)	0.5 (0.05, 7)	0.5 (0.05, 7)	0.5
540	1,080	1,600	7,000	2500
2.7	2.7	2.7	2.8	3.1
760 g	1.3 kg	1.9 kg	16 kg 20 kg wet 23 kg filled	4.9 kg 6.0 kg wet

* See section "5 Binding Capacity", page 19

4 Materials

Membrane materials

Matrix	Stabilized reinforced cellulose
--------	---------------------------------

Membrane thickness membrane volume = membrane area	275 μm 1 mL = 36.4 cm^2
---	---

Nominal pore size	> 3 μm
-------------------	-------------------

Ion exchanger ligand STIC PA	Weak anion STIC PA; primary amine (R-NH_2)
------------------------------	--

Capsule materials

Outer cage, inner core, end caps, capsule housing, nonwoven, fleece	Polypropylene
--	---------------

O-ring in vent valve (except nano)	EPDM (ethylene propylene diene monomer)
------------------------------------	--

Cassette materials

Outer cage, seal, nonwoven, fleece	ABS, silicone, polyethylene, stable to gamma irradiation
------------------------------------	---

5 Binding Capacity

Data are based on dynamic binding capacity measurements at 10% breakthrough using 3 layers of 5 cm² membrane discs (15 cm² total area membrane thickness of 275 μm) arranged in a holder and run at 10 mL/min.

Typical dynamic binding capacity 10%	Reference protein and buffer
1.4 mg/cm ² (50 mg/mL)	BSA (bovine serum albumin) in 20 mM Tris/HCl 150 mM NaCl, pH 7.5

6 Installation

The contents of the package is described in chapter “11 Ordering Information”, page 50. When unpacking capsules, protect the inlet and outlet connectors from damage. Never keep or place the capsule directly on the floor on the connectors. This might damage the sanitary adapters.

For unpacking of Jumbo 2.5 L, take the capsule including the styrene foam and end protectors out of the box and place it upright on the end protectors.

Move the Jumbo trolley (see chapter “11.2 Accessories”, page 52) in place. Then remove upper foam protection and transparent bag. Lift the Jumbo directly onto the trolley (inlet is up and the arrow imprinted on the banderole is pointing down). We recommend to connect the Jumbo with the trolley by the three screws delivered with the trolley. To ensure safe unpacking, the protective caps on inlet and outlet should stay until you use the unit. Store the caps when you plan to autoclave (see chapter “7.3 Autoclaving”, page 27). Remove before venting.

The capsules and cassettes should be installed in an upright position in the process flow. In this position the inlet is up. The flow is guided to an external channel passing through the membrane layers to an internal channel and to the outlet of the capsule (see Fig. 3). Install the capsule in-line with a prefilter (0.2 μm or 0.45 μm) in front of the device to prevent blockage or pressure build-up.

For using the cassettes you need an appropriate cassette holder and one Manifold Set (see chapter "11.2 Accessories", page 52). Before use you must read the Pilot Filter Holder manual, order no. 85037-547-72 or Process | Double Process Filter holder manual order no. 85037-553-19. If you plan to use a different filter holder from other manufacturers, you have to contact your nearest Sartorius office for technical directions.

Unpack the Manifold Set containing one inlet and one outlet plate. Place the “INLET” marked plate at one end of the holder. “THIS SIDE UP” is readable on top. Place the “OUTLET” marked plate at the other end of the holder, again “THIS SIDE UP” is readable from the top. The fluid channels of both plates are oriented to each other.

**⚠ The cassettes must be oriented in the lowest possible position in the holder otherwise the system may leak.
The cassettes must have the same lot number.**

Put the desired number of Sartobind cassettes between the manifolds (see Fig. 6). Provided “THIS SIDE UP” is readable the cassettes are correctly installed.

The maximum clamping force for 1–13 cassettes with Pilot and Process holders before start of use is 25 kilonewton or kN (200 bar, 2900 psig) for installed 12.6 cm² piston area. Now close the all DRAIN and VENT valves of the manifold plates by the pinch clamp manually.

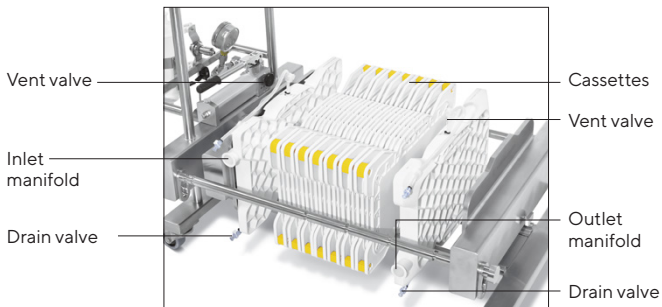


Fig. 6: Insert cassette(s) between the manifold inlet and outlet plates on the Pilot Filter Holder.

Connect the inlet and outlet plates with 1½ inch tri-clamp to the process solution. Maximum pressure for the set-up of 1 to 13 cassette(s) is 2 bar (0.2 MPa, 29 psig). Make sure that pump peak pressure caused by pulsation stays below this limit too.

7 Operation

7.1 Venting

It is important to remove air from the unit completely.

All capsules except nano carry vent valves (see Fig. 1). The vent valves are equipped with hose barb connectors for the fluid spilled out during venting. After unpacking check vent valve position. When turning anticlockwise, the valve is open, when turning clockwise, the valve is closed. Before opening the vent valve, please connect the valves with flexible tubing (inner diameter 6 mm) to waste. During venting of capsules please do not exceed 0.05 MPa (0.5 bar | 7.3 psi) pressure, as the vent valve O-ring could change its position which will result in insufficient closing of the valve.

For appropriate venting, open the vent valve screw $\frac{1}{3}$ turn to left until all air is replaced by fluid. For venting the cassettes, tubes with quick connectors are attached to the inlet and outlet manifolds and closed by a pinch clamp.

For the nano 1 mL fill a 10–20 mL Luer syringe with equilibration buffer and connect to the capsule. Hold capsule upright (outlet is up) and expel air as shown in Fig. 7. If you still detect any air in the filled unit, close the outlet, hold the syringe up and move the plunger slightly up and down that air bubbles can ascend into the syringe. Another method is to connect a second empty syringe to the top of the nano and expel air and buffer into that syringe, disconnect the upper syringe to push out air and reconnect to the nano, turn it and purge the solvent back and forth.

Very small air bubbles observed directly below the inlet of the nano do not disturb separations. The capsule will function normally as long as the small air bubbles remain outside of the membrane bed.

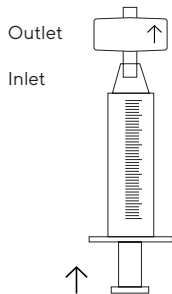


Fig. 7: Filling the Sartobind nano with a Luer syringe for air removal

7.2 Cleaning and equilibration

The devices have to be cleaned in place directly before use with 1 N NaOH, 30 min at 20°C. Preferentially work at room temperature as low temperature increases viscosity of solvents. Also cold NaOH can induce swelling of the cellulose matrix significantly.

1. For sanitization use 30 membrane volumes (MV) of 1 N NaOH solution at a flow rate of 1 MV/min.
2. Flush with 100 MV of equilibration buffer (e.g. 20 mM Tris/HCl, 150 mM NaCl, pH 7.5) at 5 MV/min.

7.3 Autoclaving

The cassette material is not compatible to autoclaving but to gamma irradiation. If you need sterile cassettes, please contact your nearest Sartorius office.

The capsules can be autoclaved once at 121°C for 30 minutes at 1 bar (0.1 MPa | 14.5 psi). Prewet the capsule with equilibration buffer. When using water you may detect a lower flow rate which is due to swelling.








The protective caps enclosed in the Jumbo delivery must be reinstalled on inlet and outlet connectors of the Jumbo. Close valves immediately after sterilisation. For autoclaving Sartobind Jumbo refer to separate autoclaving instructions enclosed in delivery.

7.4 Recommended flow rates and equilibration volumes

Membrane adsorbers can be run at much higher flow rate per volume than resin columns. The recommended flow rates for membrane adsorbers with 4 mm bed height are between 10 to 30 membrane volumes per minute. This recommendation is only a guideline since buffers and samples have different compositions and viscosities. Please test your planned flow rates using a small scale device to ensure that they fit into your pump capacities and device pressure limits. Lower flow rates than the recommended ones can also be used but will typically not improve binding capacity or overall performance. Cold room temperature increases buffer viscosity and possibly back pressure.

The equilibration volume is typically 100 membrane volumes depending on the type of buffer.

For the cassettes, flow rate and equilibration volumes have to be multiplied with the number of cassettes in use.

							
Membrane volume (MV)	1 mL	75 mL	200 mL	400 mL	600 mL	2.5 L	800 mL
Rec. flow rate (L/min)	0.02	1.5	4	8	12	50	16**
Rec. equilibration volume* (L)	0.1	7.5	20	40	60	250	80**

* Refer to 7.2 Cleaning and equilibration

** Multiply with number of used cassettes

7.5 Buffer conditions

Ionic strength of buffers during loading can be much higher than for conventional anion exchange membrane adsorbers up to 20 mS/cm. The pKa of the chosen buffer should not exceed ± 0.5 pH units of the operation pH. It should be filtered with 0.2 μm or 0.45 μm filters before use and the quality of water and chemicals should be of high purity.

⚠ It is recommended to use monovalent buffers e.g. TRIS or Acetate. Multivalent buffers like phosphate or citrate will reduce binding capacity for proteins. On the other hand, contaminants such as DNA and endotoxin can still bind depending on character of multivalent buffer up to ~150 mM to achieve even better separation of target molecules from contaminants. Application of pure water may lead to a reversible swelling of the membrane and may reduce permeability.

7.6 Selection of pH and salt conditions

In ion exchange chromatography a charged molecule is bound to oppositely charged groups attached to the insoluble matrix. This binding is reversible by application of salt ions to the buffer eluting the molecule. The pH value at which a biomolecule has no net charge is the isoelectric point: pI . If the pH of the buffer is below the isoelectric point (rule of the thumb at least 1 pH unit) a protein for example carries a positive net charge and will bind to a cation exchanger (Sartobind S). If the pH of the buffer is above its iso-electric point (at least 1 pH unit), it will bind to anion exchangers (Sartobind Q or Sartobind STIC PA).

Consider the character of the amine ligand of Sartobind STIC PA as a weak anion exchanger. It means that approaching higher pH the positive charge lowers. To optimize the binding capacity and load volume, multiple pH values should be tested (e.g. on 96 well plates).

Conventional ion exchangers are loaded at low conductivity. Proteins are easily eluted by adding e.g. 1 M NaCl. Sartobind STIC PA binding is also influenced by salt but higher levels of salt are needed to elute the molecules. At a level of e.g. 150 mM NaCl where conventional anion exchanges do not bind, the salt tolerant membrane shows good binding capacity. To remove protein from the membrane, higher salt concentration than for conventional ion exchanger shall be needed.

7.7 Contaminant removal from therapeutic proteins and other sources in flow-through mode

For contaminant removal from products such as monoclonal antibodies, pH conditions in the range of pH 6 to 8 should be used. Contaminants include highly negatively charged DNA, endotoxins, protein contaminants, some host cell proteins and viruses. The product of interest, the monoclonal antibody with isoelectric points (pI) of 8 – 9.5 for example, will not bind and pass through the Sartobind STIC PA. The influence of the flow rate is very low.

7.8 Sample preparation

The sample should be adjusted to the starting buffer conditions and be prefiltered through a 0.2 µm membrane filter e.g. Sartopore® 2 capsule. For small volumes in the mL range use a 0.2 µm Minisart® filter with Luer outlet (order number 16532-K for polyethersulfone or 16534-K for cellulose acetate membrane).

⚠ Unfiltered feed will block the Membrane Adsorber and lead to capacity loss and increased back pressure. We recommend inline filtering during operation. With increase of pressure replace filter and restart.

7.9 Washing

When using capsules in bind & elute mode, wash with equilibration buffer after sample loading.

7.10 Elution

To elute the target protein use buffer with appropriate salt concentration. Take into account that STIC has been developed for single-use and that the original binding capacity typically cannot be restored due to strong binding of the primary amine ligand to the negatively charged species.

7.11 Draining

You may drain the capsules by application of air or nitrogen pressure (<1 bar| 14.5 psi) to the inlet of the device.

⚠ A dual air regulator system is recommended to prevent over-pressurization of the Sartobind device. The first regulator should reduce line air pressure to 2 bar. The second regulator, positioned immediately upstream of the Sartobind capsule, should reduce the 2 bar regulated supply pressure to the <1 bar (14.5 psi) for a capsule and 0.5 bar (7.3 psi) for 1 up to 13 cassettes draining pressure.

7.12 Regeneration and Storage

After elution and wash with equilibration buffer a regeneration step with 1 N NaOH for 1 hour may clean the STIC but consider as described in the preceding chapter, that binding capacity typically cannot be restored completely. If necessary store in 20% ethanol in equilibration buffer.

7.13 Chemical stability

The devices are stable against all commonly used buffers in chromatography. Do not use oxidizing agents.

7.14 Operation of the Sartobind nano with peristaltic pumps or liquid chromatography (LC) systems

After the unit is filled completely with equilibration buffer, close the outlet of the Sartobind nano and remove the syringe. Start the LC system or peristaltic pump at a low flow rate. When fluid emerges, stop the pump, connect the tubing to the inlet of the Sartobind nano. Make sure that no air is introduced. Remove the cap from outlet.

Run the pump until fluid emerges from the outlet of the unit and stop it. Then connect the outlet of the unit via Luer adapter to the LC detector and proceed with loading. If your system pressure is too high, refer to your LC system manual to remove any flow restrictor after the UV cell, as the system may generate a pressure above the allowed maximum pressure. As membrane adsorbers run typically at much higher flow rates than columns, there is no risk of bubble formation in the UV cell when removing the restrictor.

7.15 Scaling up

Complete the break through experiments for the target compound to be bound on the membrane matrix. After optimisation of binding conditions for the contaminants, the purification step can be scaled up to a larger capsule.

Recommendations:

Maintain

- Bed height (automatically kept constant when using capsules with same bed height)
- Linear flow (automatically kept constant when using capsules with same bed height)
- Sample concentration

Increase (see scaling factors in the following table)

- Sample loading volumes
- Volumetric flow rate
- Membrane volume

Scale up is done preferably by keeping the bed height and changing the the membrane volume as the calculation is most simple. Other methods for scale up via residence time will lead to same result. The residence time is the membrane volume divided by the flow rate.


Using the Sartobind nano 1 mL the scale up factor for flow rate and binding capacity is equal to a multiplication factor of the membrane volumes of the listed scale up devices:

Size	Membrane volume [mL]	Factor to increase* (from nano)
nano	1 mL	-
5"	75 mL	75
10"	200 mL	200
20"	400 mL	400
30"	600 mL	600
Jumbo	2.5 L	2500
Cassette	800 mL	800
Cassettes**	10.4 L	10,400

* Flow rate and binding capacity;

** 13 Cassettes as example

Example: After breakthrough experiments with the nano you found 500 fold binding capacity is needed. You will choose the 600 mL capsule. Then increase flow rate by a factor of 600 mL/min too.

 **Keep sample concentration constant in lab and production scale. Watch out for volumes in the piping and flow rate in whole system.**

8 Integrity Test by Diffusion

The integrity of the membrane adsorber can be controlled by a diffusion test.

The testing procedure describes the diffusion test for pre and post use. The test is intended to discriminate between defective and intact devices and to detect major bypasses, large holes and faulty assembly.

8.1 Installation for test

Test procedure has been generated with current Sartocheck® instrument family e.g. Sartocheck® 4 Plus (26288), 4 (16288) or 3 Plus (16290).


Use of earlier Sartocheck® instruments will generate faulty data. Install capsule as shown in Fig. 8.

Please note that the test procedure with other vendor's integrity testers can require a different set up.

8.2 Operation of test

8.2.1 Pre-washing of device

Pre-wash with 30 membrane volumes (MV) of buffer or 0.9% NaCl in water at recommended flow rate.

-  The capsule needs to be pre-washed with the testing solvent, to remove any glycerol. The washing solution should be at room temperature. Keep the unit in an upright position for proper venting and open the vent screw on top of the device until all air is replaced by testing solvent.

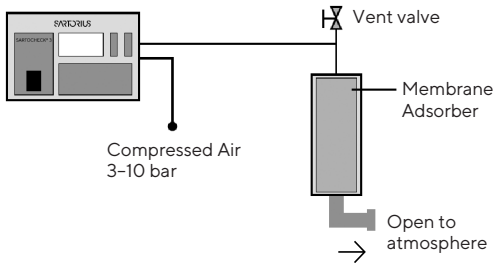


Fig. 8: Set up of diffusion test with Sartochek® 4 Plus, 4 or 3 Plus

8.2.2 Diffusion measurement with Sartochek® 4 Plus

- Choose **Programming** in the main menu
- Choose **Diffusion Test**

Choose the test pressure, stabilization and testing time for your device from the table. If you set the Net Volume to zero, Sartochek® automatically measures the upstream void volume including tubing.

Test parameters

Size	Bed height (mm)	Membrane volume (MV)	Test pressure mbar (psi)	Stabilisation time (min)	Testing time (min)	Diffusion max. mL/min
nano	4 mm	1 mL	100 (1.45)	2	1	-
5"	4 mm	75 mL	100 (1.45)	2	1	75
10"	4 mm	200 mL	100 (1.45)	3	1	75
20"	4 mm	400 mL	100 (1.45)	3	1	75
30"	4 mm	600 mL	100 (1.45)	3	1	75
Jumbo	4 mm	2.5 L	100 (1.45)	3	1	150
Cassette(s)*	4 mm	0.8-10.4 L	100 (1.45)	5	1	15-195

* Diffusion max. per 4 mm cassette is 15 mL/min multiplied by number of cassettes

8.2.3 Results and evaluation for one device

- Diffusion \leq Diffusion max.:
Test passed (diffusion value on the print out)
- Diffusion $>$ Diffusion max.:
Test failed (red text on the print out)

9 Troubleshooting

Problem	Possible cause	Action
Air bubbles can be seen	Incomplete air removal	Small air bubbles seen in the top of the unit do not interfere with the purification as long as they do not touch the membrane bed. If too much air is enclosed, repeat removal as described in chapter "7.1 Venting", page 24.
I installed the capsule upside down	Installation of capsule may be easier in the process flow	Validation has been done with a process flow from top to bottom. Thus it is clearly recommended to use capsules in the described flow direction (Feed enters capsule on top and leaves it on bottom).
I deviated from the CIP and flushing equilibration procedure		The capsules have been qualified and validated according the given procedure. If a deviation is necessary, the results may also deviate from the given validation data.
High back pressure	Material has not been filtered	Prefilter with 0.2 µm or 0.45 µm filter before processing through the unit (preferentially inline).

Problem	Possible cause	Action
High back pressure	Material has been filtered but was stored before purification	Proteins can form aggregates within hours or during operation. Thus we recommend to prefilter inline by attaching a 0.2 µm filter in front of the adsorber. When you observe again pressure built up, replace the filter.
	LC system generates high pressure	Remove restrictor after the UV cell.
	The adsorber is clogged membrane fouling	Replace unit. You may backflush within given flow and pressure limits, perform a regeneration cycle.
	Viscosity swelling effects	Work at room temperature, avoid lower temperatures
	Pure water leads to swelling of membrane	Add sodium chloride or use ionic buffers

Problem	Possible cause	Action
I have to use water for wetting the capsules	Economic considerations; easier usage	You may use pure water but you have to expect a decreased flow rate which may lead to a higher backpressure. The use of water will not change diffusion testing values. The flow rate returns to normal after using buffer again.
Target molecule is not bound	Conditions for binding are insufficient	Process conditions, e.g. prefiltration, pH, conductivity, multivalent buffers etc. have to be checked and optimized. Sartobind STIC membrane is salt tolerant.
Binding capacity is not sufficient	Process conditions not optimized	Use larger adsorber device, or: connect two adsorbers (same size) in series (i.e connect outlet of first adsorber to inlet of second) to achieve higher binding capacity. As a rule of thumb the pressure doubles when the flow rate is kept constant and the number of membrane layers is doubled.
Incomplete elution	Strong binding	Use capsule only for polishing in flowthrough mode.

Problem	Possible cause	Action
I want to reuse the STIC capsules	Economic reasons	The ligand strongly binds the contaminants and cleaning with 1 N NaOH does typically not restore 100 % of the binding capacity depending on the character of your sample and the contaminants. Sartobind STIC has been developed for mainly single-use application to avoid the revalidation.
I need to remove DNA endotoxins but I have to use PBS	I heard PBS is a multivalent buffer which reduces binding capacity	Go ahead with the PBS as DNA is a high negatively charged species and will bind to STIC. The same accounts for endotoxins. You may even work at pH conditions which normally would not be accessible as target protein could be bound.
A vertical line is seen on one capsule side when filled	Membrane edge visible	No action necessary. It can be visible the edge of the fleece touching the inner tube.
I purged with air or nitrogen and lost flow and binding capability.	Air has entered into the pores	See troubleshooting “Applied bubble point instead of diffusion test” below.

Problem	Possible cause	Action
Accidentally a bubble point test instead of diffusion test has been run	Operation error	The membrane has then to be purged extensively to remove all the air which has been pressed into the pores. If properly purged, the diffusion test can be run successfully and the device works as expected.

10 Quality Assurance

This product is tested for protein dynamic binding capacity and flow rate. Sartobind membranes have been tested for protein dynamic binding capacity, flow rate, thickness, and evenness. Capsules, cassettes and membranes are manufactured in a controlled environment. The product meets all Sartorius Stedim Biotech standards for traceability, production and specifications as given here or exceeded them as certified in the quality assurance certificate enclosed. A validation and an extractables guide are available on request.

11 Ordering Information

11.1 Products

Order number	Description and type of connectors	Quantity
96STPA42DN-11-A	Sartobind STIC PA nano 1 mL, 4 mm, Luer female connectors, 2 PEEK adapters Luer male to UNF 10-32 female, manual, certificate	4
96STPA42D9MFF--A	Sartobind STIC PA 75 mL, 4 mm, ¾" sanitary clamp, manual, certificate	4
96STPA42D1GSS	Sartobind STIC PA 200 mL, 4 mm, 1½" sanitary clamp, manual, certificate	1
96STPA42D2HSS	Sartobind STIC PA 400 mL, 4 mm 1½" sanitary clamp, manual, certificate	1
96STPA42D3KSS	Sartobind STIC PA 600 mL, 4 mm, 1½" sanitary clamp, manual, certificate	1




Order number	Description and type of connectors	Quantity
96STPA42D3NSS	Sartobind STIC PA Jumbo 2.5 L, 4 mm 1½" sanitary clamp, 2 protective caps, manual, autoclaving instructions, certificate	1
98STPA42D-L	Sartobind STIC PA Cassette 0.8 L, 4 mm, 1½" sanitary clamp via manifold set (accessory), manual, certificate	1

11.2 Accessories

Order number	Description	Quantity
1ZA---0004	Adapter Luer male to UNF 10-32 female, PEEK	1
1ZAOGV0003	Adapter UNF 10-32 female to sanitary ¾", 25 mm, polyoxymethylene	2
5ZGI--0001	Holder for 1 × 200 to 1,200 mL (10-30") capsule, stainless steel, 3 legs	1
5ZALB-0002	Distribution adapter for 3 × 200 (10-30") to 1200 mL capsules, 1 × 2", 3 × 1½", sanitary, stainless steel	1
7ZAL-V0013	Reducing adapter 1½" (50.5 mm) to ¾" (25 mm), sanitary	1
7ZAL-V0010	Reducing adapter 2" (64 mm) to 1½" (50.5 mm), sanitary	1
9ZGL--0102	Trolley for Jumbo 2.5 or 5 L, stainless steel	1
16290	Sartocheck 3 Plus Integrity Tester	1
26288---FT	Sartocheck 4 Plus Filter Integrity Tester	1

Order number	Description	Quantity
29Z-S00001	Manifold set for Sartoclear® Sartobind®, 1½" sanitary clamp	2
2ZGL--0005	Pilot filter holder for Sartoclear® Sartobind®	1
2ZGL--0006	Process filter holder for Sartoclear® Sartobind®	1
2ZGL--0007	Double process filter holder for Sartoclear® Sartobind®	1
2ZGL--0008	Drip pan for Pilot Filter holder	1
2ZGL--0015	Drip pan for Process and double Process Filter Holder	1

12 Dimensions and Connections

Membrane volume 4 mm bed height	 1 mL	 75 mL	 200 mL
Size	nano	5"	10"
Dimensions in mm	37×33 H×Ø	190×77 H×Ø	350×100 H×Ø
Connectors	Luer female	Sanitary ¾" 25 mm outer, 14 mm inner diameter"	Sanitary 1½" 50.5 mm outer, 36 mm inner diameter
Gaskets	n.a.	¾", inner diameter 16 mm	1½", inner diameter 35.8 mm
Recommended internal diameter of flexible tube for hose barb version	n.a.	12.7 mm	n.a.

n.a.=not available



400 mL

20"

570×100
H×Ø

Sanitary 1½"
50.5 mm outer,
36 mm inner
diameter

1½",
inner diameter
35.8 mm

n.a.



600 mL

30"

810×100
H×Ø

Sanitary 1½"
50.5 mm outer,
36 mm inner
diameter

1½",
inner diameter
35.8 mm

n.a.



2.5 L

Jumbo

850×302
H×Ø

Sanitary 1½"
50.5 mm outer,
36 mm inner
diameter

1½",
inner diameter
35.8 mm

n.a.



0.8 L

Cassette

634×387×49
W×L×Ø

Via manifold:
Sanitary 1½"
50.5 mm outer,
36 mm inner
diameter

For manifold:
1½", inner diameter
35.8 mm

n.a.

Sartorius Stedim Biotech GmbH
August-Spindler-Strasse 11
37079 Goettingen, Germany

Phone: +49 551 308 0
www.sartorius.com

First published:
December 15, 2016
Sartorius Stedim Biotech GmbH,
Goettingen, Germany

Printed in the EU on paper bleached
without chlorine.
DIR: 2624862-000-00

The information and figures
contained in these instructions
correspond to the version date
specified below.

Sartorius reserves the right to make
changes to the technology, features,
specifications and design of the
equipment without notice.

Masculine or feminine forms
are used to facilitate legibility in
these instructions and always
simultaneously denote the other
gender as well.

Copyright notice:

This instruction manual, including
all of its components, is protected
by copyright. Any use beyond the
limits of the copyright law is not
permitted without our approval.
This applies in particular to
reprinting, translation and editing
irrespective of the type of media
used.

© Sartorius Germany

Last updated:
08 | 2020